

Description: HighTaq DNA Polymerase is a novel genetically engineered version of *Taq* DNA polymerase which provides high robustness against inhibitors (e.g. bio-molecules like polysaccharides, fats and proteins). The performance of the enzyme ensures high sensitivity and reliable amplification. The polymerase replicates DNA at 72 °C. It catalyses the polymerization of nucleotides into duplex DNA in 5' → 3' direction and possess a 5' → 3' exonuclease activity. High Taq DNA Polymerase is available as high concentrated solution with 75 U/μl.

Content

Ref No.	S111105 (5 U/μl)	111105 (5 U/μl)	111125 (5 U/μl)	111106 (75 U/μl)	color
HighTaq DNA Polymerase	Sample size	500 units	2500 units	1000 units	blue
Incomplete NH ₄ * Reaction Buffer (10x)	1.8 mL	1.8 mL	5x 1.8 mL	2x 1.8 mL	red
Complete NH ₄ ** Reaction Buffer (10x)	1.8 mL	1.8 mL	5x 1.8 mL	2x 1.8 mL	yellow
Complete KCl*** Reaction Buffer (10x)	1.8 mL	1.8 mL	5x 1.8 mL	2x 1.8 mL	black
MgCl ₂ 100 mM	1 mL	1 mL	5x 1 mL	2x 1 mL	green
Datasheet	1	1	1	1	--

* Incomplete NH₄ Reaction Buffer (10x): 160 mM (NH₄)₂SO₄, 670 mM Tris-HCl (pH 8,8), 0.1 % Tween 20

** Complete NH₄ Reaction Buffer (10x): 160 mM (NH₄)₂SO₄, 670 mM Tris-HCl (pH 8,8), 0.1 % Tween 20, 25 mM MgCl₂.

*** Complete KCl Reaction Buffer (10x): 500 mM KCl, 100 mM Tris-HCl (pH 8,8), 0.1 % Tween 20, 15 mM MgCl₂.

Applications: HighTaq DNA Polymerase is suitable for all regular applications but especially for those where inhibitors might decrease the performance. HighTaq DNA Polymerase effectively directs PCR with the template up to 5 kb in length. High Taq DNA Polymerase high concentrated (75 U/μl) is recommended for lyophilization.

Concentration: 5 units/μL

Sensitivity: detection of ≥ 6 DNA molecules

Unit definition: One unit of activity is defined as the amount of enzyme required to incorporate 10 nmoles of dNTP into acid-insoluble DNA fraction in 30 minutes at 72°C.

Additionally provided: 1 tube MgCl₂ (100 mM)

Recommended MgCl₂ concentration: 1.5 mM – 6 mM

Quality control

- 98% protein homogeneity in 10% SDS-PAGE
- No detectable exo-/endonuclease activities
- PCR amplification tests with different templates
- PCR amplification tests without templates as Negative Control
- “no template” test with primers complementary to a region auf 16S bacterial ribosomal genes

Storage condition: -20 °C

Pipetting scheme and thermocycler protocol:

Components	Volume / 50 μ L PCR-Reaction	Final concentration
10 x PCR-Buffer	5 μ L	1x
dNTP-Mix (40 mM)	1 μ L	800 μ M (200 μ M each)
Upstream Primer	variable	0.1 - 0.5 μ M
Downstream Primer	variable	0.1 - 0.5 μ M
HighTaq DNA Polymerase	0.25 - 1.0 μ L	1.25 - 5.0 units
Template DNA	variable	10 to 500 ng/reaction
Sterile dest. water	Adjust to 50 μ L final volume	

Separate $MgCl_2$ solution can be used for optimization. If incomplete buffer is used **titrate $MgCl_2$** for optimal PCR results with following recommendation (see table):

Final $MgCl_2$ conc. [mM]	1.5	2	2.5	3	3.5	4	4.5	5	5.5	6
Volume [μ L] of 100 mM $MgCl_2$ / 50 μ L	0.75	1	1.25	1.5	1.75	2	2.25	2.5	2.75	3

Thermocycler protocol

step	time	temperature
initial denaturation	2 minutes	94 °C
30 cycles:		
denaturation	10 seconds	94 °C
annealing	20 seconds	55 - 68 °C *
extension	1 minute	72 °C

* Usually the optimal annealing temperature is 5 °C below the melting temperature of the primers.

Notes:

Program the cycler according to the manufacturer's instructions.

Each program should start with an initial denaturation step at 94 °C for 2 to max. 5 min.

Recommended elongation time: 1 min/ 1 kb of target.

For maximum yield and specificity, temperatures (annealing) and cycling times should be optimized for each new template target or primer pair.

We Bring The World Of Biotechnology To You



www.bocascientific.com

(781) 686-1631

info@bocascientific.com